


# PRODUCT EXPLODED VIEW

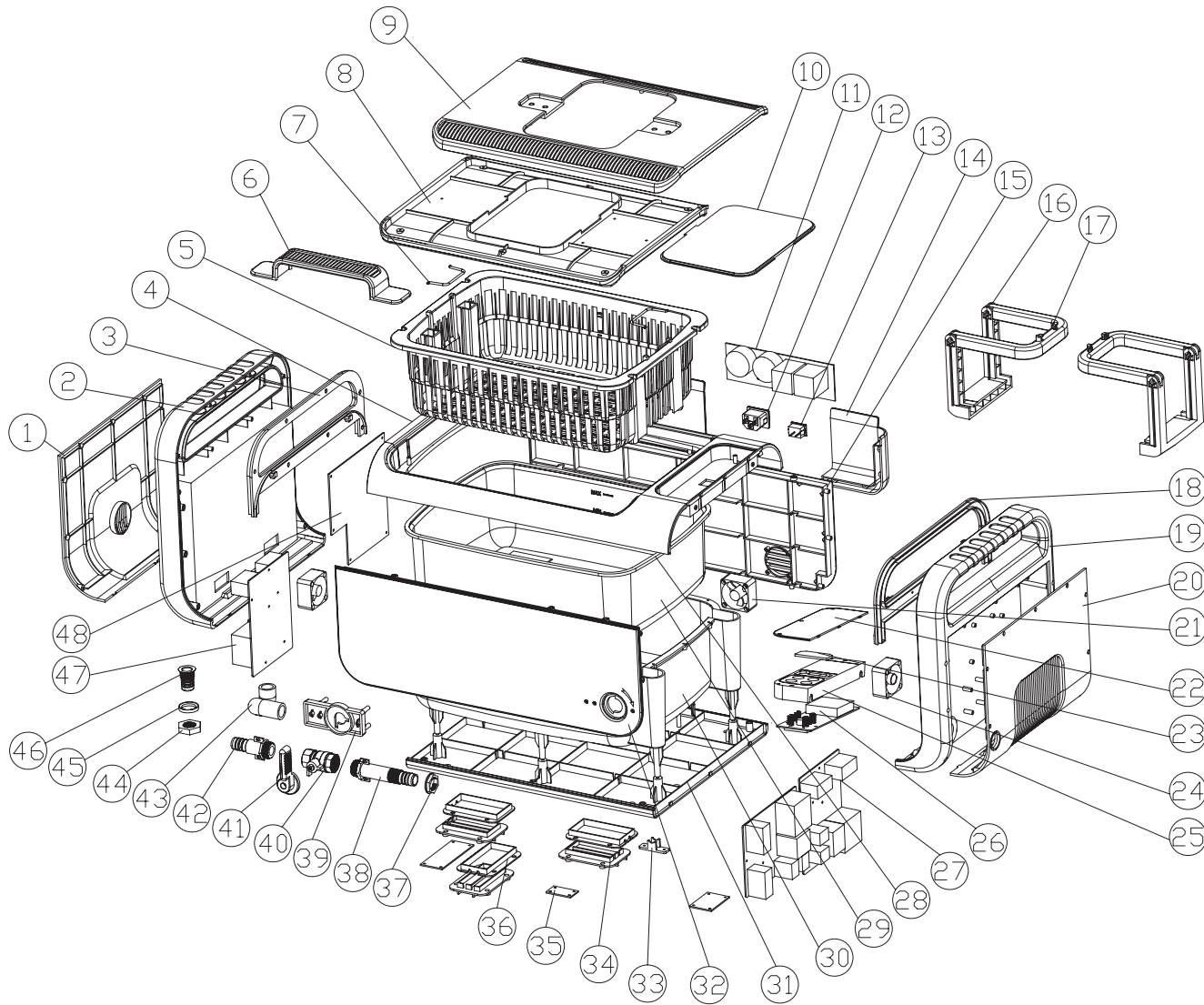

No.	Item	Material
①	Left cover	ABS
②	Left panel/handle	ABS
③	Left inside cover	ABS
④	Top ring	ABS
⑤	Plastic basket	ABS
⑥	Top handle	ABS
⑦	Stainless steel handle	SUS 304
⑧	Lid-bottom	ABS
⑨	Lid-top	ABS
⑩	Lid-window	AS
⑪	Power switch PCB	94V <sub>0</sub>
⑫	Power socket	Standard part
⑬	Power switch	Standard part
⑭	Lid holder	ABS
⑮	Back cover	ABS
⑯	Basket foldable handle 1	ABS
⑰	Basket foldable handle 2	ABS
⑱	Right inside cover	ABS
⑲	Right panel/handle	ABS
⑳	Right cover	ABS
㉑	Cooling fan	Standard part
㉒	Control panel cover	PMMA
㉓	Illumination window	AS
㉔	Cooling fan	Standard part
㉕	Control panel box	AS
㉖	LED	Fire retardant ABS
㉗	Control PCB	94V <sub>0</sub>
㉘	Silicon seal	Silicone rubber
㉙	Stainless steel tank	SUS 304
㉚	Inner tank	ABS
㉛	Bottom cover	ABS
㉜	Front cover	ABS
㉝	Thermocouple holder	ABS
㉞	Heater cover	PA 66
㉟	Insulation	94V <sub>0</sub>
㊱	Heater pocket	PA 66
㊲	Nut	ABS
㊳	Spout	ABS
㊴	Bracket	ABS
㊵	Ball valve	Standard part
㊶	Valve handle	ABS
㊷	Elbow	ABS
㊸	Hose	Silicone rubber
㊹	Nut	AL
㊺	Washer	AL
㊻	Spout	SUS 304
㊼	Transducer PCB	94V <sub>0</sub>
㊽	Power PCB	94V <sub>0</sub>